Vangerowstraße 14
69115 Heidelberg
+49 (0)6221 - 90 50 50

www.finescience.de
europe@finescience.de

DEUTSCH	(ab S. 2)
ENGLISH	(from p. 3)
FRANÇAIS	(à partir de la p. 4)
ITALIANO	(da p. 5)
ESPAÑOL	(desde p. 6)

FST # 18000-45 & 18000-50

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INSTRUCTIONS

SURGICAL INSTRUMENT STERILIZER FST No. 18000-45 AND -50

Fill instrument with glass beads provided, to within 3 mm below rim.

Insert electrical plug into wall socket.

Unit will take 15 to 20 minutes to reach operating temperature of 250°C.

Carefully insert the working part of clean surgical instrument into well after unit has reached operation temperature and leave for 10 to 20 seconds. Do not leave in longer as instrument becomes too hot to handle. Sterilization in on contact with beads, so tubes (canulae) cannot effectively be sterilized.

Unit can be left in the "ON" position all day, and is safe for use in laminar flow cabinets.

Only when unit is switched off and cool, cover with lid to reduce dust accumulation. Do not cover with lid when unit is hot.

Should unit overheat and trip safety switch, let it cool off completely, then turn it over and push red button visible inside the unit as viewed through one of the cooling holes.

The glass beads should be cleaned routinely and replaced in a dry state.

Use a spoon for removal of the glass beads, since otherwise there is a slight chance of beads getting stuck in the rim of the well.

IMPORTANT: This unit should only be used for research purposes. DO NOT use in situations which involve humans or valuable animals. DO NOT USE in surroundings where inflammable gases/substances are involved.

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