
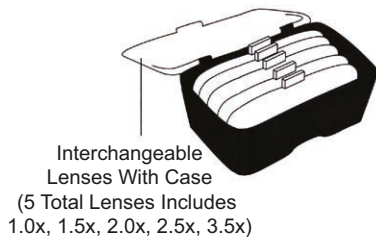


**Features & Contents**  
**Item No. 28080-10**


**Interchangeable Magnifier Headset**

# Features & Contents

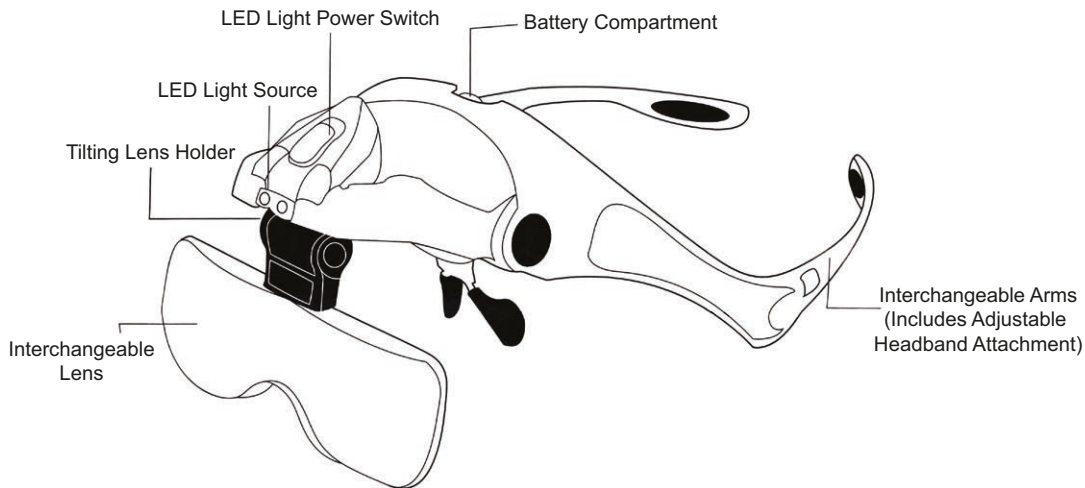

### Changing the Batteries

1. Remove the back panel by pushing the tab.

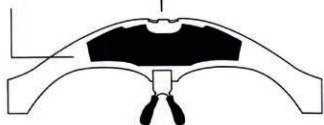


2. Insert (3) AAA batteries and replace back panel.\*




*\*Batteries Not Included*

### Changing Lenses

1. Choose the desired lens and push into the bottom of the tilting lens holder.

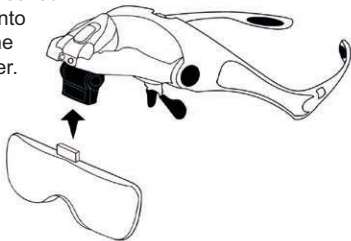


2. To remove the lens, pull the lens from the bottom of the tilting lens holder.

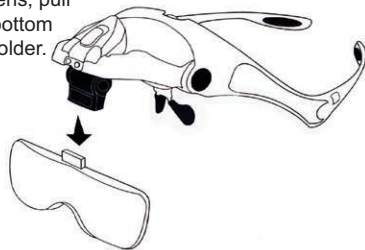


### Replacing the Arms with the Headband

1. Push the tab located at the start of the arm. Pull the arm out while pushing the tab in.

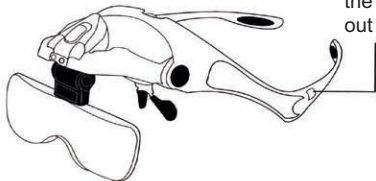

2. Insert the end of the headband into the tab. The tab will click when fully inserted. Repeat step one to remove the headband.

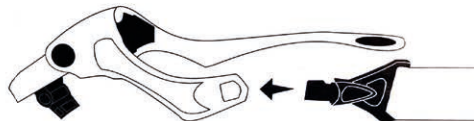

**Fine Science Tools (USA) Inc.**

4000 East 3rd Avenue, Suite 100  
Foster City, CA 94404-4824  
USA

**Telephone** +1 800-521-2109 / +1 650-349-1636

**Fax** +1 650-349-3729

**E-Mail** [info@finescience.com](mailto:info@finescience.com)

**Web** [finescience.com](http://finescience.com)

**Fine Science Tools (CAN) Inc.**

410-197 Forester Street  
North Vancouver, British Columbia  
Canada V7H 0A6

**Telephone** +1 800-665-5355 / +1 604-980-2481

**Fax** +1 800-665-4544 / +1 604-987-3299

**E-Mail** [canada@finescience.com](mailto:canada@finescience.com)

**Web** [finescience.ca](http://finescience.ca)

**Fine Science Tools GmbH**

Vangerowstraße 14  
69115 Heidelberg  
Germany

**Telephone** +49 62 21 - 90 50 50

**Fax** +49 62 21 - 90 50 590

**E-Mail** [europa@finescience.com](mailto:europa@finescience.com)

**Web** [finescience.de](http://finescience.de)