# Instruction Manual for Demagnetizer

Item No. 18000-15

EN

DE

FR

ES

IT

PT

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JA

KO

This laboratory accessory is perfect for demagnetizing forceps, screwdrivers, as well as other microsurgical instruments and parts. It can also magnetize tools, such as forceps, which can come in handy when working with small screws that can disappear when dropped. This magnetizer has a high power to weight ratio and is housed in a non-breakable plastic case. Due to the high power of this unit, the current should not be kept on continuously, but in case it is left on, this unit features an automatic shut-off.

### **Contents:**

- 1 Demagnetizer with line cord
  - 120 V A.C., 60Hz., 300 watts

### **NOTE:**

Please read these instructions carefully and completely before operating the Demagnetizer. Ensure unit is properly plugged into a grounded electrical outlet of proper voltage and amperage, and ensure no liquids come in contact with the unit at any time. Accident or injury may occur if these precautions are not followed.

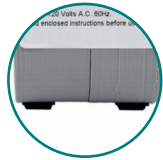
# PRODUCT FEATURES:

One Push Button Operation


8 Feet Cord Length

Rubber Feet to stabilize unit

## TO DEMAGNETIZE:

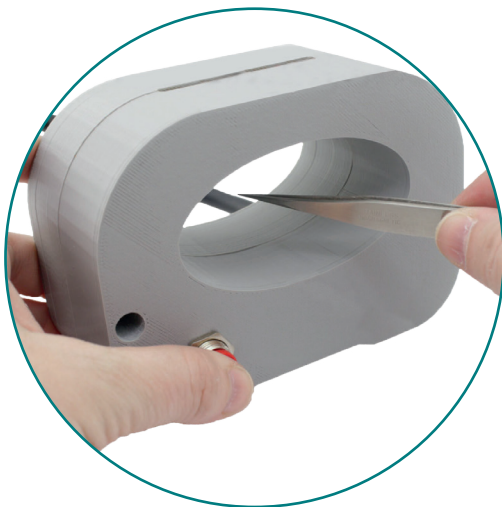- 1.** Hold the item to be demagnetized inside the center of the housing opening. Use caution to ensure the part is not touching the housing.
- 2.** Press and hold the button and withdraw the item slowly in a straight line from the housing opening for approximately three inches, keeping the item in line with the center of the housing opening, then release the button.
- 3.** The item should now be demagnetized. Repeat procedure as necessary.

## TO MAGNETIZE:

- 1.** Position the item to be magnetized inside and touching the sides of the housing opening.
- 2.** Press and hold the button for four or five seconds then release the button.
- 3.** The item should now be magnetized. Repeat procedure as necessary.

## DUTY CYCLE:

- This unit has superior field strength due to the relatively high amperage it draws. The consequence is a fairly rapid heating of the field coil. It is recommended that a single duty cycle be limited as follow: 5 seconds on and 3 minutes off, or three 5 second cycles on then 10 minutes off.
- This unit is wired with a thermal protector. If the above duty cycle are not observed, the thermostat may open and the unit will not operate until it has cooled off sufficiently. This unit is not intended for continuous duty but for intermittent use only.

## STERILIZATION:

- GAD, EtO only
- Do not steam autoclave

## CONTACT INFO:

### **Fine Science Tools (USA) Inc.**

4000 East 3rd Avenue, Suite 100  
Foster City, CA 94404-4824

#### **USA**

Telephone: +1 800-521-2109 / +1 650-349-1636

Fax: +1 800-523-2109/+1 650-349-3729

E-Mail: [info@finescience.com](mailto:info@finescience.com)

Web: [finescience.com](http://finescience.com)

### **Fine Science Tools (CAN) Inc.**

410-197 Forester Street  
North Vancouver, British Columbia

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### **Fine Science Tools GmbH**